

# Ringbio




## Ringbio Milk Rapid Test Kit

Accurate

Sensitive

Reliable

internationally approved

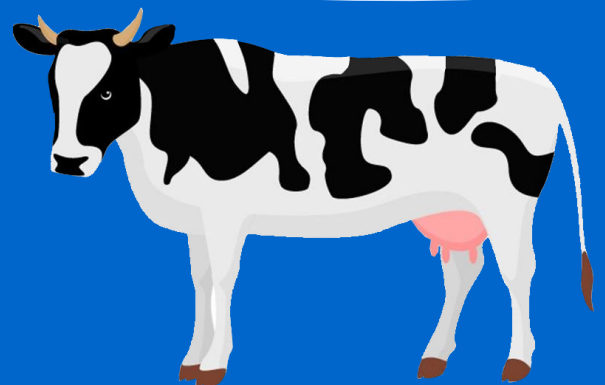

*Accurate, Sensitive, Reliable and Internationally approved tests*



These tests can be used to monitor antibiotic residues and mycotoxins, melamines, etc in milk and various food samples. Ringbio is also working with ILVO and other international laboratories to validate the tests.

Product Code	Product	Approval
100002	$\beta$ -lactams & tetracyclines, BT Combo Test Kit	ILVO, Russia, Czeck, Turkey, Iran, etc
100015	$\beta$ -lactams & tetracyclines & sulfa drugs, BTS TriTest 3in1	Argentina, ILVO in preparation
100029	$\beta$ -lactams & tetracyclines & sulfa & quinolones, BTSQ QuaTest 4in1	ILVO
100018	$\beta$ -lactams & tetracyclines & streptomycin & CAP, BTSC QuaTest 4in1	Russia, ILVO in preparation
100025	Aminoglycosides, GNKS QuaTest 4in1	ILVO, Russia
100024	Quinolones, macrolides, lincomycin, erythromycin, QMLE QuaTest 4in1	Russia
100004	Aflatoxin M1 Rapid Test Kit	ILVO
420003	Camel milk fraud test kit	ILVO
R6002	Folic Acid Rapid Test Kit	ILVO
E6001	Bovine Lactoferrin ELISA Kit	ILVO in preparation



## About ILVO

ILVO is short for Flanders Research Institute for Agriculture, Fisheries and Food in Belgium, which is a world leading research insititute on food safety. It is the AOAC and AFNOR expert laboratory for chemical analysis.

## *$\beta$ -lactams & tetracyclines, BT Combo Test Kit*



A lateral flow rapid test kit based on receptor assay to detect multiple  $\beta$ -lactams & tetracyclines antibiotics in 10min.

This kit is validated by



ILVO



Russian VNIMI  
Institute of Dairy Industry



Scan this code to view the detailed information on [ringbio.com](http://ringbio.com)

### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

### *Order information*

Product Code	Reaction Model	Detection Level
100002	5'+5', room temp.	EU MRL
102002	5'+5', room temp.	EU MRL
104002	3'+7', room temp.	Russian MRL
105002	3'+7', room temp.	EU MRL
106002	3'+3', 40°C	EU MRL

## *$\beta$ -lactams & tetracyclines & sulfa drugs, BTS TriTest 3in1*



A lateral flow rapid test kit based on receptor assay to detect multiple  $\beta$ -lactams & tetracyclines & sulfa antibiotics in 6min.

This kit is validated by



Argentina INTI

### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

### *Order information*

Product Code	Reaction Model	Detection Level
100015	3'+3', room temp.	EU MRL
101015	3'+5', room temp.	EU MRL
102015	3'+5', room temp.	Beta EU MRL, TC Russian MRL
103015	3'+3', room temp.	EU MRL, higher sensitivity

ILVO validation in preparation.



Scan this code to view the detailed information on [ringbio.com](http://ringbio.com)

## *$\beta$ -lactams & tetracyclines & sulfa & quinolones, BTSQ QuaTest 4in1*



A lateral flow rapid test kit based on receptor assay to detect multiple  $\beta$ -lactams & tetracyclines & sulfa antibiotics in 6min.

This kit is validated by



### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

### *Order information*

Product Code	Reaction Model	Detection Level
100029	3'+7', room temp.	EU MRL
101029	3'+7', room temp.	TC Russian MRL, other EU MRL



Scan this code to view the detailed information on [ringbio.com](http://ringbio.com)

## *$\beta$ -lactams & tetracyclines & streptomycin & CAP, BTSC QuaTest 4in1*



A lateral flow rapid test kit based on receptor assay to detect multiple beta-lactams, tetracyclines, streptomycin & chloramphenicol antibiotics in 10min.

This kit is validated by



Russian VNIIM  
Institute of Dairy Industry

ILVO validation in preparation



Scan this code to view the detailed information on [ringbio.com](http://ringbio.com)

### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

### *Order information*

Product Code	Reaction Model	Detection Level
100018	3'+7', room temp.	EU MRL
101018	3'+7', room temp.	Ultra sensitive LOD
102018	3'+7', room temp.	TC EU MRL

## *Aminoglycosides, GNKS QuaTest 4in1*



A lateral flow rapid test kit based on receptor assay to detect multiple gentamicin, neomycin, kanamycin & streptomycin antibiotics in 10min.

This kit is validated by



ILVO



Russian VНИИ  
Institute of Dairy Industry



Scan this code to view the detailed information on [ringbio.com](http://ringbio.com)

### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.

### *Order information*

Product Code	Reaction Model	Detection Level
100025	3'+7', room temp.	EU MRL
101025	3'+7', room temp.	Ultra sensitive LOD

## *Quinolones, macrolides, lincomycin, erythromycin, QMLE QuaTest 4in1*



A lateral flow rapid test kit based on receptor assay to detect multiple quinolones, macrolides, lincomycin, erythromycin antibiotics in 10min.

This kit is validated by



Russian VNIMI  
Institute of Dairy Industry

### *Kit Component*

- 12 tubes of rapid tests, 8 strips per tube.
- Microwell holder
- Plastic pipette or micropipette tips,
- Kit instruction

### *Samples*

- Raw milk, UHT milk, pasteurized milk, etc.
- Water, meat, egg, honey, etc.


### *Order information*

Product Code	Reaction Model	Detection Level
100024	3'+7', room temp.	EU MRL
101024	3'+7', room temp.	Ultra sensitive LOD



Scan this code to view the  
detailed information on  
[ringbio.com](http://ringbio.com)



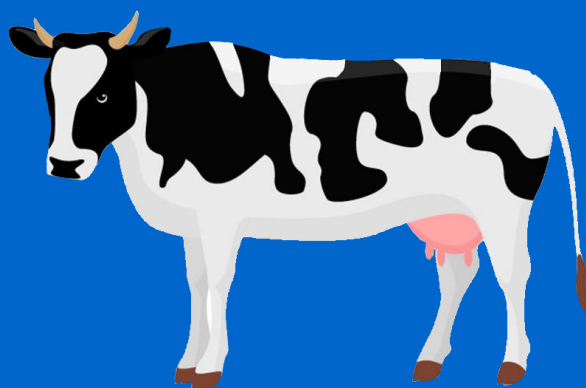

# Ringbio Milk Rapid Test Kit

**Ring Biotechnology Co Ltd**

**Add:** Building 3, Zhongtongtai Science Park,  
No. 11, Kechuang 14th St,  
Beijing 100176, CHINA

**Phone:** +86.10.56267496

**Web:** [www.ringbio.com](http://www.ringbio.com)

**Email:** [info@ringbio.com](mailto:info@ringbio.com)